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# User Guide

[www.wtavet.com.br](http://www.wtavet.com.br)

**Aquecedores de Tubos – WTA**

**Tubes Warmers - WTA**

## ADVERTÊNCIA

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## Index

<b>WARNING</b> .....	<b>13</b>
Copyright.....	13
Comments .....	13
<b>PRESENTATION</b> .....	<b>14</b>
<b>ACCESSORIES</b> .....	<b>14</b>
<b>APPLICATION</b> .....	<b>15</b>
<b>TECHNICAL FEATURES</b> .....	<b>15</b>
<b>FUNCTION SPECIFICATION</b> .....	<b>16</b>
<b>IMPORTANT PRECAUTIONS</b> .....	<b>17</b>
Repairs: .....	17
Cares and Transport: .....	17
Cleaning: .....	17
Falling Objects Inside the Equipment:.....	17
Power Cable:.....	17
Rain and Thunderstorms: .....	17
<b>INSTALATION, SETTING AND OPERATION</b> .....	<b>18</b>
Connecting the Equipment .....	18
Setting the Temperature .....	18
<b>WARRANTY</b> .....	<b>18</b>
Voiding the Warranty.....	19
<b>PROBLEM SOLUTION</b> .....	<b>20</b>
Malfunction in the Electrical Network .....	20
Heating the Appliance .....	21
<b>TECHNICAL ASSISTANCE SERVICE / CUSTOMER SERVICE</b> .....	<b>21</b>

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## USER GUIDE

### WARNING

To prevent accidents and/or damage, read and follow carefully this instruction guide before using the product.

Warning for dangerous voltage presence.

Warning for important operating instructions.

### Copyright

Tubes Warmers - WTA user guide

WTA, Cravinhos, State of São Paulo, Brazil

The content of this guide is protected by copyright laws and cannot be reproduced, distributed, disclosed neither stored without prior written permission of WTA - Watanabe Tecnologia Aplicada Eireli - EPP.

In order to obtain authorization, additional information or to clarify doubts, please contact our Technical Assistance Department by e-mail, letter or phone.

Phone: +1(979) 324-6168

e-mail: [usa1@wtatech.com](mailto:usa1@wtatech.com)

\*Opening hours: From Monday to Friday, 8:30 am to 5:30 pm (GMT -03:00)

©WTA - Watanabe Tecnologia Aplicada Eireli - EPP - All rights reserved.

### Comments

The information, illustrations and specifications contained in this user guide assume the latest features of the product available at the time of its publication.

Due to the use of new technologies and raw materials, WTA reserves the right to change the design and/or technical specifications of the product always as necessary and without prior notice, without incurring in any obligations.

This manual is an integral part of the product. Therefore, the consumer should save it for any queries.

## PRESENTATION

Congratulations on your choice. You have just purchased a WTA product, company which adds quality and technology in what it manufactures.


WTA is a Brazilian company specialized in solutions for animal reproduction. Our products are developed for working with cattle, sheep and horses, with recognized approval by the most renowned specialists in the world.

Structured in anodized aluminum offer an elegant design and, conditions that facilitate both internal and external cleaning.

For best product performance we advise you to read this manual.

## ACCESSORIES

When purchasing your equipment, make sure your accessories are intact and included in the package:

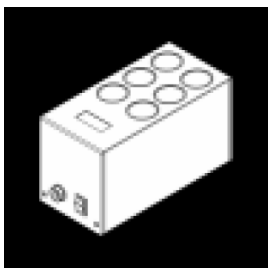Tubes Warmer

Power Cable

Power Supply

**Note:** If any accessory is missing or damaged, contact the WTA customer service department within 3 (three) days of the product delivery date.

## APPLICATION

Tubes Warmers are equipment that provides thermal safety for procedures that require heating of biological material and laboratory solutions.

Practical and easy to use, the equipment has an electronic sensor that maintains the temperature while working with  $\pm 1^{\circ}\text{F}$  ( $0.5^{\circ}\text{C}$ ) tolerance of variance.

## TECHNICAL FEATURES

- Manufactured in aluminum.

- Capacity:

  - AT50-i: 6 tubes of 50mL.

  - ATM: 4 tubes of 50 mL + 4 tubes of 15 mL + 6 microtubes.

  - 1(50) / 10(15): 1 tube of 50 mL + 10 tubes of 15 mL.

  - 15(15): 15 tubes of 15 mL.

- Approximate weight: 7,45 lb (3,400kg).

- Dimensions (W x H x H): 200 x 92 x 112mm.

- Maximum power: 64.5W.

- AC power source: 100~240Vac - 1.4A - 50/60Hz.


- DC unit supply: 15 Vdc, 4.3A.

- Temperature range: 68.0 to 113.0°F (20.0 to 45.0°C).

- Temperature variation:  $\pm 1^{\circ}\text{F}$  ( $0.5^{\circ}\text{C}$ ).

- Product warranty: 3 years.

## FUNCTION SPECIFICATION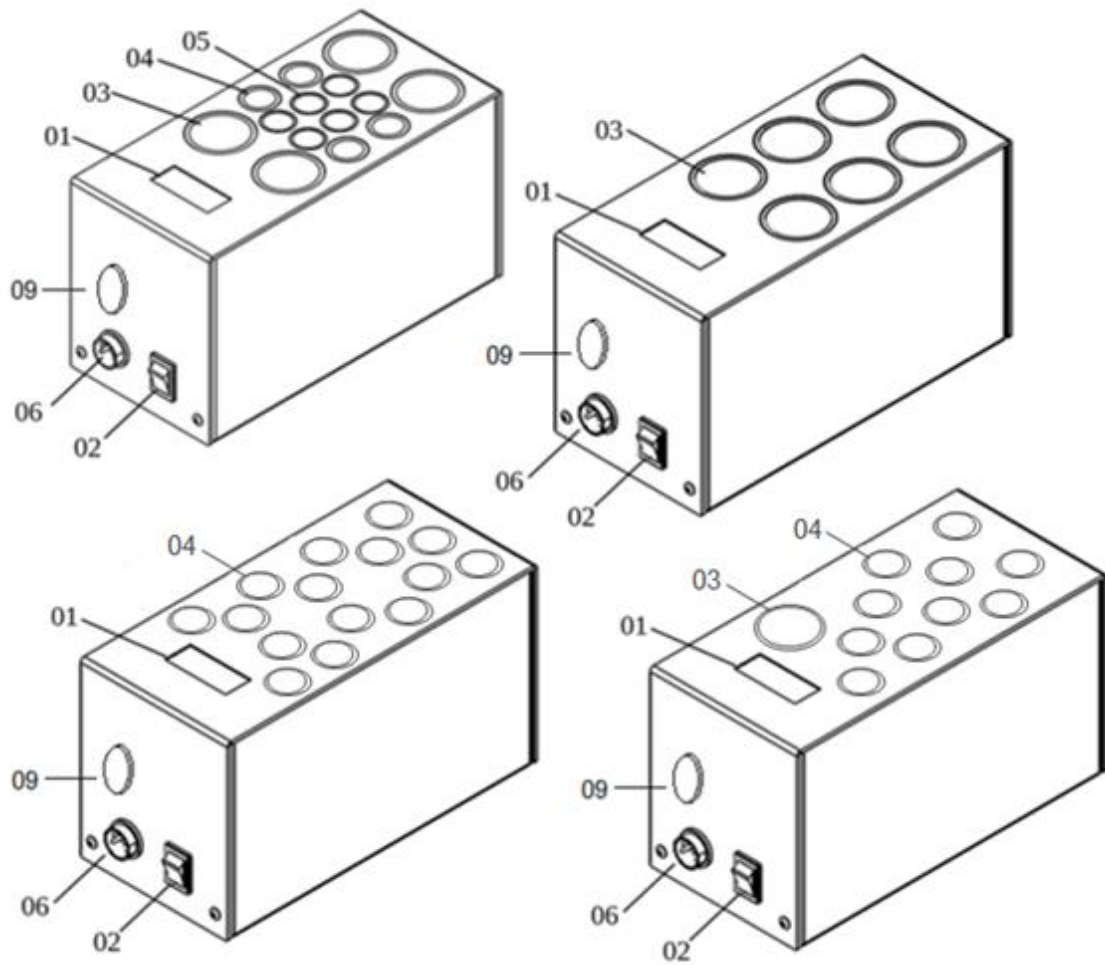

01 - LCD display

02 - Settings button

03 - 50 mL tube warmer

04 - 15 mL tube warmer


05 - Cryotube warmer

06 - Power input DC

07 - AC power supply adapter

08 - DC output

09 - On/off button

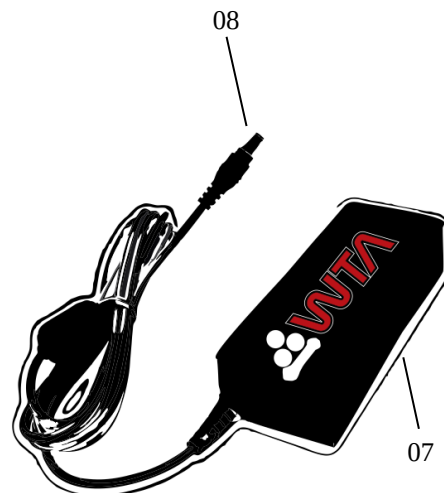## IMPORTANT PRECAUTIONS

### Repairs:

Do not open the device. All repairs and maintenance should be performed by authorized professionals or by WTA service.

### Cares and Transport:

Avoid shaking during transport which can damage it. Make sure that the equipment is properly seated and protected against any accidents.

### Cleaning:

Clean the exterior with a soft cloth slightly moistened with water. Do not use alcohol, thinner or other solvents as they will cause damage device's finishing.

### Falling Objects Inside the Equipment:

Do not place recipients or cups containing liquid, or coins, paper clips or other metallic objects on the control module. If any liquid or any objects falls inside the equipment, turn off the power immediately and contact the service center.

### Power Cable:

Do not step on the power cable or put any kind of objects (furnishing, etc.) over it.

### Rain and Thunderstorms:

In storms with electric discharge events disconnect the equipment from the AC power.

## INSTALATION, SETTING AND OPERATION

### Connecting the Equipment

For safety insert the power supply connector (08) into the equipment input (06). Then connect the power supply (07) to the power network 100 ~ 240Vac using the power cable.

### Setting the Temperature

When the power supply (07) is connected to the power network (and/or press on/off button (09)) a beep sounds and the display (01) will show the equipment name, serial number and model. Then another beep will be emitted and the display will show SETPOINT. Press the "+" or "-" (02) buttons to adjust the desired temperature. The set value will be stored in memory for future use.

The temperature can be adjusted to a range of 68.0 to 113.0°F (20.0 to 45.0°C).

This equipment will not provide cooling if temperature setted was lower than ambient temperature.

To turn off press on/off button (09) (or remove the plug from the wall outlet).

## WARRANTY

WTA - Watanabe Applied Technology Eireli - EPP follows strict internal quality control standards in the products manufacture. In addition to the legal guarantee of 90 (ninety) days, provided for in the Consumer Protection Code, it guarantees its customers an additional 33-month Warranty, counted from the date of issue of the purchase invoice, provided that the equipment has been installed and according to the guidelines contained in the User's Manual.

Even if ownership of the equipment is transferred, the warranty will continue to be valid according to the terms and conditions specified above.

This product has been carefully evaluated at every stage of the manufacturing process. This warranty will always accompany you, valid only in the country where the product was purchased.

It is understood by Additional Warranty, the free repair of the product and / or the replacement of parts that, according to technical opinion of the WTA, present a manufacturing defect.

Only WTA Technical Assistance professionals are entitled to repair defects for the guarantee, by presenting the sales tax (original) to the first buyer / consumer.

### **The Warranty Service**

- Assistance in WTA technical assistance.
- Parts used for maintenance, subject to the exceptions provided in the item "Exclusion of Warranty".

The product should be sent directly to the WTA observing that:

- The product must be properly packed, if possible in its original packaging, together with the manual, cables and other accessories that have been delivered at the time of purchase.
- The product will be pre-evaluated by WTA technicians prior to repair.
- Shipping, freight and packaging costs will only be charged to WTA during the legal guarantee period (90 days). After this period such costs will be the customer's responsibility.

### **Voiding the Warranty**

This warranty does not cover defects or defects arising from the following situations:

- Installation and use of the product in an improper way or in disagreement with the respective guidelines and information described in this Manual, including installation, external adjustment, cleaning and use of accessories not in accordance with technical specifications.
- Defects caused by accidents, falls, atmospheric phenomena (rains, floods, electric discharges), excessive humidity (water, liquids in general), salinity, sinister (theft or theft), loss or improper transport of the product.
- Installation, adjustment, repair or servicing by persons outside WTA service.
- Invoice and / or serial number adulterated or withdrawn.

- Modifications, alterations or violation of the original product.
- Product connected to the mains improper or subject to excessive fluctuations.
- The product is purchased in balance, showcase or demonstration, this being characterized by means of identification in the invoice of purchase and / or price lower than the one practiced in the market.

**Important:** WTA is not liable for damage to equipment during transportation. When purchasing your product make sure on the delivery act that it is not defective or with apparent defects. If you notice any irregularities, return the product immediately.

### **This Warranty Statement Does Not Include**

- Periodic reviews.
- Displacement and expense of technician by authorized WTA service.

### **Responsibility**

The WTA has no control or influence over the conditions under which the equipment is used, as well as the methodologies employed. Thus, WTA assumes no responsibility for the results, use and / or performance of the equipment (including: damages for loss of profits, interruption of business activity, loss of business information or other pecuniary losses). WTA expects that the equipment will be used by trained personnel. The purchaser acknowledges that the equipment is subject to all terms and conditions described in this manual.

## **PROBLEM SOLUTION**

Occasionally many apparent defects may be something simple to solve. If you notice any problems with the equipment, check the following procedures.

### **Malfunction in the Electrical Network**

- Check that the power cable is connected correctly.
- Disconnect it from the mains, wait 5 seconds and connect again.

## **Heating the Appliance**

- Check that the temperature is set correctly.
- Check that the cables are properly connected.
- Disconnect them from the mains, wait 5 seconds and connect again.

If the problem persists, contact your service representative.

## **TECHNICAL ASSISTANCE SERVICE / CUSTOMER SERVICE**

If there is any doubt or if it is necessary additional information about warranty, assembly, installation, adjustment or utilization of the product, please contact our Customer Service Department:

Opening hours: From Monday to Friday, 8:30 am to 5:30 pm (GMT -03:00)

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Images contained in the Guide are merely illustrative.

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Todos os direitos reservados | All rights reserved